



## ZYCEINE STERILE SPORICIDE FLUID

HYPOCHLOROUS ACID

ITEM NO 9420 (200ML X 15 PER CASE)

EFFECTIVE CLEANROOM SPORICIDES ARE USUALLY AGGRESSIVE CHEMICALS AND WHILST THEY CAN GIVE GOOD MICROBIAL EFFICACY THIS CAN OFTEN LEAD TO CORROSION PROBLEMS AND DAMAGE TO EXPENSIVE EQUIPMENT.

ZYCEINE IS A UNIQUE TWO COMPONENT SPORICIDE THAT WHEN ACTIVATED GIVES GOOD MICROBIAL EFFICACY. ZYCEINE IS A TOTALLY STABLE PRODUCT UNTIL ACTIVATED.

ZYCEINE HAS A PROVEN FUNGICIDAL EFFICACY OF GREATER THAN A 3-LOG REDUCTION IN 2 MINUTES, AND A YEASTICIDAL EFFICACY GREATER THAN A 4-LOG REDUCTION IN 2 MINUTES.

- FAST ACTING SPORICIDE.
- EFFICACY INDEPENDENTLY VALIDATED.
- WFI QUALITY WATER – HIGH PURITY WATER.
- 0.2 MICRON FILTERED.
- FULLY TRACEABLE WITH SUPPORTING ANALYSIS.
- GAMMA IRRADIATION FOR STERILITY.
- 2 YEAR UNOPENED SHELF LIFE.
- ALSO AVAILABLE IN 1L FLUID AND 68GSM LOW PARTICULATE POLYCELLULOSE WIPES



### DIRECTION FOR USE:

AGMA ZYCEINE IS A UNIQUE SPORICIDE, MANUFACTURED TO HELP CLEANROOM OPERATORS MEET THE STANDARDS OF ANNEX 1. TO ACTIVATE, ROLL THE POUCH FROM THE BOTTOM TOWARDS THE FRANGIBLE SEAL. AS PRESSURE BUILDS, SQUEEZE GENTLY UNTIL THE MID-POINT SEAL BREAKS. LEAVE TO REST FOR 5 MINUTES TO FULLY ACTIVATE BEFORE USING.

Use Biocides Safely. Always read the label and product information.

THE INFORMATION CONTAINED IN THIS DATA SHEET IS ACCURATE TO THE BEST OF OUR KNOWLEDGE AND BELIEF. NO WARRANTY IS, HOWEVER, GIVEN OR IMPLIED. CONSULT AGMA LTD CONCERNING ALL NON-STANDARD APPLICATIONS.

ISSUE NO. 1-03/25