



# ZYCEINE STERILE SPORICIDE WIPES (DB)

HYPOCHLOROUS ACID

ITEM NO. 0131 (10 X 5 WIPES)

EFFECTIVE CLEANROOM SPORICIDES ARE USUALLY AGGRESSIVE CHEMICALS AND WHILST THEY CAN GIVE GOOD MICROBIAL EFFICACY THIS CAN OFTEN LEAD TO CORROSION PROBLEMS AND DAMAGE TO EXPENSIVE EQUIPMENT.

ZYCEINE IS A UNIQUE TWO COMPONENT SPORICIDE THAT WHEN ACTIVATED GIVES GOOD MICROBIAL EFFICACY. ZYCEINE IS A TOTALLY STABLE PRODUCT UNTIL ACTIVATED.

ZYCEINE HAS A PROVEN FUNGICIDAL EFFICACY OF GREATER THAN A 3-LOG REDUCTION IN 2 MINUTES, AND A YEASTICIDAL EFFICACY GREATER THAN A 4-LOG REDUCTION IN 2 MINUTES.

- FAST ACTING SPORICIDE.
- EFFICACY INDEPENDENTLY VALIDATED.
- MADE WITH WFI QUALITY WATER.
- 0.2 MICRON FILTERED.
- MANUFACTURED IN GRADE C/D CLEANROOMS.
- RE-SEALABLE TO PREVENT DRYING OUT.
- 68GSM LOW PARTICULATE POLYCELLULOSE WIPE.
- 28CM X 38CM FOR LARGE SURFACE AREA COVERAGE.
- BROAD SPECTRUM ACTIVITY.
- FULLY TRACEABILITY WITH SUPPORTING ANALYSIS.
- GAMMA IRRADIATED FOR STERILITY.
- 2 YEAR UNOPENED SHELF LIFE.
- DOUBLE BAGGED.
- ALSO AVAILABLE IN 200ML AND 1L FLUID POUCHES.



## DIRECTION FOR USE:

AGMA ZYCEINE IS A UNIQUE SPORICIDE, MANUFACTURED TO HELP CLEANROOM OPERATORS MEET THE STANDARDS OF ANNEX 1. TO ACTIVATE, ROLL THE POUCH FROM THE BOTTOM TOWARDS THE FRANGIBLE SEAL. AS PRESSURE BUILDS, SQUEEZE GENTLY UNTIL THE MID-POINT SEAL BREAKS TO TRANSFER THE FLUID ONTO THE WIPES. AGITATE THE POUCH THOROUGHLY FOR 30 SECONDS, LEAVE TO REST FOR 30 SECONDS – REPEAT OVER A 5 MINUTE INTERVAL.

Use Biocides Safely. Always read the label and product information.

THE INFORMATION CONTAINED IN THIS DATA SHEET IS ACCURATE TO THE BEST OF OUR KNOWLEDGE AND BELIEF. NO WARRANTY IS, HOWEVER, GIVEN OR IMPLIED. CONSULT AGMA LTD CONCERNING ALL NON-STANDARD APPLICATIONS

ISSUE NO. 1-01/26