



STERILE ALKALINE DETERGENT WIPES (DB)

IN WFI PRE-IMPREGNATED – READY TO USE
ITEM 0125 (10 X 10 WIPES)

FOR EFFECTIVE DISINFECTION TO BE ACHIEVED IT IS IMPERATIVE THAT SURFACES ARE FREE FROM ANY VISIBLE SOILING OR BUILD-UP OF RESIDUE FROM PREVIOUS DISINFECTANT APPLICATIONS.

AGMA HAS CONSIDERABLE EXPERIENCE IN SURFACTANT TECHNOLOGY AND STERILE CLEANING / DISINFECTION. WE HAVE COMBINED THIS EXPERTISE TO DEVELOP STERILE ALKALINE DETERGENT WIPES, SPECIFICALLY FOR USE IN CLEANROOMS AND PHARMACIES.

WITH HIGHLY EFFECTIVE SURFACE CLEANING PROPERTIES, AGMA STERILE ALKALINE DETERGENT IS IDEAL WHEN A DEEP CLEAN IS REQUIRED.

- pH12 IDEAL FOR DEEP CLEANS AND HIGH PROTEIN LOAD CLEANING.
- PARTICULARLY EFFECTIVE AGAINST PROTEINACEOUS AND ORGANIC RESIDUES.
- REMOVES BUILD-UP THAT NEUTRAL OR ACIDIC CLEANERS MAY NOT.
- LOW FOAMING.
- GAMMA IRRADIATED FOR STERILITY.
- 0.2 MICRON FILTERED.
- MADE WITH WFI WATER.
- MANUFACTURED IN GRADE C/D CLEANROOMS.
- QMS TO BS EN ISO 9001:2015.
- MANUFACTURED IN ACCORDANCE WITH THE cGMP GUIDELINES.
- DOUBLE BAGGED.
- FULLY TRACEABLE WITH SUPPORTING ANALYSIS.



DIRECTION FOR USE

AGMA ALKALINE DETERGENT WIPES ARE MANUFACTURED TO HELP CLEANROOM OPERATORS MEET THE STANDARDS OF ANNEX 1. IDEAL FOR USE ON CLEANROOM HARD SURFACES TO REMOVE RESIDUES AND FOR USE BETWEEN ROTATING OF DISINFECTANTS. AN INITIAL DEEP CLEAN SHOULD BE PERFORMED PRIOR TO THE USE OF ANY DISINFECTANT TO HELP THE REMOVAL OF ANY SOILS OR HEAVY RESIDUE BUILD UP. USE AGMA'S HIGH PURITY WATER TO WIPE DOWN AFTER USE.

THE INFORMATION CONTAINED IN THIS DATA SHEET IS ACCURATE TO THE BEST OF OUR KNOWLEDGE AND BELIEF. NO WARRANTY IS, HOWEVER, GIVEN OR IMPLIED. CONSULT AGMA LTD CONCERNING ALL NON-STANDARD APPLICATIONS.

ISSUE NO. 1-01/26