



STERILE POLYESTER DRY WIPES (DB)

140GSM POLYESTER DRY WIPES – READY TO USE ITEM 1280 (10 X 25 WIPES)

STERILE POLYESTER DRY WIPES HAVE VERY LOW LEVELS OF PARTICLES AND FIBRES, SO THEY REDUCE CONTAMINATION RISK. THEIR SMOOTH, DURABLE SURFACE MAKES THEM IDEAL FOR WIPING DOWN SENSITIVE EQUIPMENT AND SURFACES WITHOUT SNAGGING OR LEAVING FIBRES.

PAIRING OUR 100% POLYESTER STERILE DRY WIPES WITH AGMA'S RANGE OF CLEANROOM COMPATIBLE STERILE FLUID DISINFECTANTS ALLOWS THE REMOVAL OF MICROBIAL CONTAMINANTS AND INHIBITS THEIR GROWTH.

STERILE POLYESTER WIPES HAVE THE ADDED BENEFIT OF BEING ABLE TO PHYSICALLY REMOVE PARTICLES AND SPORES WHILE PROVIDING GOOD DISINFECTION AGAINST VEGETATIVE MICROBES.

- 100% HIGH-QUALITY POLYESTER WIPE.
- LASER SEALED EDGE FOR REDUCED FIBRE CONTAMINATION.
- FOR USE IN GRADE A/B CLEANROOMS.
- GAMMA IRRADIATED FOR STERILITY.
- 23CM X 23CM - 140GSM WIPES.
- MANUFACTURED IN GRADE C/D CLEANROOMS.
- FULLY TRACEABLE WITH SUPPORTING ANALYSIS.
- DOUBLE-BAGGED (DB) PLUS BOX LINER.
- 25 WIPES PER PACK.
- 2-YEAR SHELF LIFE.



DIRECTION FOR USE:

AGMA STERILE POLYESTER DRY WIPES HELP CLEANROOM OPERATORS MEET THE STANDARDS OF ANNEX 1, FOR USE ON CLEANROOM EQUIPMENT AND HARD SURFACES.

APPLY THE DISINFECTANT/CLEANER (E.G., IPA) DIRECTLY ONTO THE WIPE, NOT THE SURFACE, TO PREVENT AEROSOLISING PARTICLES. THE WIPE SHOULD BE DAMP, NOT DRIPPING.

WIPE IN ONE DIRECTION USING PARALLEL, OVERLAPPING STROKES TO ENSURE YOU DON'T MISS ANY SURFACE AREAS. START AT THE CLEANEST AREA AND MOVE TO THE DIRTIEST.

Use Biocides Safely. Always read the label and product information.

THE INFORMATION CONTAINED IN THIS DATA SHEET IS ACCURATE TO THE BEST OF OUR KNOWLEDGE AND BELIEF. NO WARRANTY IS, HOWEVER, GIVEN OR IMPLIED. CONSULT AGMA LTD. CONCERNING ALL NON-STANDARD APPLICATIONS.

ISSUE NO. 1-04/26