



# STERILE NEUTRAL DETERGENT (DB)

140GSM PRE-IMPREGNATED POLYESTER WIPES – READY TO USE  
ITEM 0144 (7 X 25 WIPES)

A 100% KNITTED POLYESTER WIPE, IMPREGNATED WITH STERILE NEUTRAL DETERGENT WITH WATER FOR INJECTION (WFI) QUALITY WATER.

POLYESTER CLEANROOM WIPES HAVE VERY LOW LEVELS OF PARTICLES AND FIBRES, SO THEY LOOK TO MITIGATE CONTAMINATION RISK. THEIR SMOOTH, DURABLE SURFACE MAKES THEM IDEAL FOR WIPING DOWN SENSITIVE EQUIPMENT AND SURFACES WITHOUT SNAGGING OR LEAVING FIBRES.

WITH HIGHLY EFFECTIVE SURFACE CLEANING PROPERTIES, AGMA STERILE NEUTRAL DETERGENT IS IDEAL WHEN A DEEP CLEAN IS REQUIRED. FORMULATED TO BE LOW FOAMING IT DOES NOT LEAVE SIGNIFICANT RESIDUES OR SMEARS. IN ADDITION, ITS PH HELPS PROTECT MORE DELICATE WORK SURFACES AND THOSE PRONE TO CORROSION.

- 100% HIGH-QUALITY POLYESTER WIPE.
- LASER-CUT AND SEALED EDGE FOR REDUCED FIBRE CONTAMINATION.
- FOR USE IN GRADE A/B CLEANROOMS.
- GAMMA IRRADIATED FOR STERILITY.
- 0.2 MICRON FILTERED.
- 23CM X 23CM - 140GSM WIPES.
- RE-SEALABLE TO PREVENT DRYING OUT.
- WFI QUALITY WATER – HIGH PURITY WATER.
- MANUFACTURED IN GRADE C/D CLEANROOMS.
- FULLY TRACEABLE WITH SUPPORTING ANALYSIS.
- DOUBLE-BAGGED (DB) PLUS BOX LINER.
- 2-YEAR SHELF LIFE.



## DIRECTION FOR USE

AGMA NEUTRAL DETERGENT WIPES ARE MANUFACTURED TO HELP CLEANROOM OPERATORS MEET THE STANDARDS OF ANNEX 1. IDEAL FOR USE ON CLEANROOM HARD SURFACES TO REMOVE RESIDUES AND FOR USE BETWEEN ROTATING OF DISINFECTANTS. AN INITIAL DEEP CLEAN SHOULD BE PERFORMED PRIOR TO THE USE OF ANY DISINFECTANT TO HELP THE REMOVAL OF ANY SOILS OR HEAVY RESIDUE BUILD UP.

THE INFORMATION CONTAINED IN THIS DATA SHEET IS ACCURATE TO THE BEST OF OUR KNOWLEDGE AND BELIEF. NO WARRANTY IS, HOWEVER, GIVEN OR IMPLIED. CONSULT AGMA LTD CONCERNING ALL NON-STANDARD APPLICATIONS.

ISSUE NO. 1-06/26